

References:

1. Heymann PW, Chapman MD, Aalberse RC, Fox JW, Platts-Mills TAE. Antigenic and structural analysis of group II allergens (Der p II and Der f II) from house dust mites (*Dermatophagoides* spp.). *J Allergy Clin Immunol* 1989;83:1055-1067.
2. Lombardero M, Heymann PW, Platts-Mills TAE, Fox JW, Chapman MD. Conformational stability of B cell epitopes on Group I and Group II *Dermatophagoides* spp. allergens. *J Immunol* 1990;144:1353-1360.
3. DeBlay F, Heymann PW, Chapman MD, Platts-Mills TAE. Airborne dust mite allergens: comparison of Group II allergens with Group I mite allergen and cat allergen, Fel d I. *J Allergy Clin Immunol* 1991;88:919-926.
4. Ovsyannikova IG, Vailes LD, Li Y, Heymann PW, Chapman MD. Monoclonal antibodies to Group II *Dermatophagoides* spp. allergens: Murine immune response, epitope analysis and development of a two-site ELISA. *J Allergy Clin Immunol* 1994;94:537-546.
5. Platts-Mills TAE, Vervloet D, Thomas WR, Aalberse RC, Chapman MD. Indoor Allergens and Asthma: Report of the Third International Workshop. *J Allergy Clin Immunol*, 1997; 100:S1



www.inbio.com

Indoor Biotechnologies, Inc.
700 Harris Street,
Charlottesville, VA 22903
United States

Tel: (434) 984-2304
Fax: (434) 984-2709
Email: mail@inbio.com

Indoor Biotechnologies Ltd.
Vision Court
Caxton Place
Cardiff, Wales CF23 8HA
United Kingdom

Tel: +44 (0) 29 2167 4640
Email: Info@indoorbiotech.co.uk

Indoor Biotechnologies,
India Private Limited
Bangalore Bioinnovation Centre,
BioTech Park, Electronic City Phase 1
Bangalore-560100, India

Tel: +91-9901722009
Email: info@inbioindia.com




Mite Group 2 ELISA kit (1D8/7A1)

Product Code: EL-D2

Lot Number: XXXXX

Sample Curve:

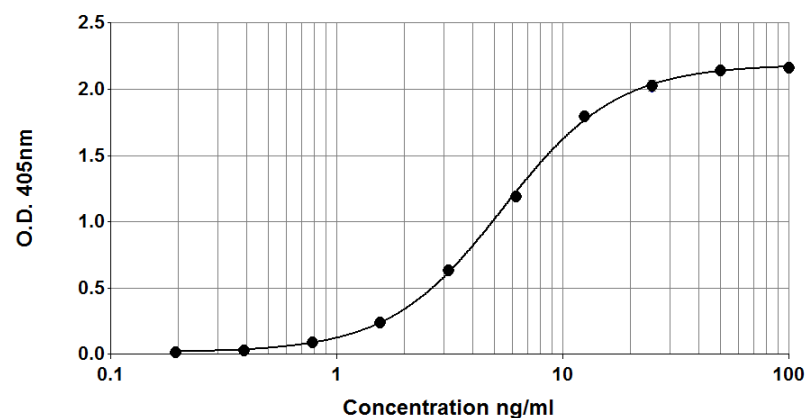

Content:

Vial 1 (red top) 100 µL
Monoclonal antibody 1D8

Vial 2 (white top) 400 µL
Universal Allergen Standard
Concentration: 1000ng/ml Der p 2

Vial 3 (brown) 100 µL
Biotinylated monoclonal antibody 7A1
Dilute: 1:1000 for use

Storage: The ELISA kit should be stored at 4°C

For research and commercial use in vitro: not for human in vivo or therapeutic use.

An InBio® product

Certificate of Analysis

Monoclonal Antibody: 1D8 (clone 1D8 C4 D8)
Immunogen: Der p 2
Isotype: Mouse IgG2A
Specificity: Binds to an epitope on dust mite *Dermatophagoides* Group 2 allergens, Der p 2 and Der f 2.
Purification: Produced in ascites and purified by ammonium sulphate precipitation and affinity chromatography using Protein A. Single heavy and light chain bands on SDS-PAGE.
Concentration: 1.35 mg/ml in phosphate buffered saline, pH 7.4
 Based on A280 for IgG (1.42=1mg/ml), 0.22µm filtered, preservative free.
Lot Number: xxxxx

Monoclonal Antibody: Biotin 7A1 (clone 7A1 H1 G3)
Immunogen: Der f 2
Isotype: Mouse IgG1
Specificity: Binds to an epitope on dust mite *Dermatophagoides* Group 2 allergens, Der p 2 and Der f 2.
Purification: Produced in ascites and purified by affinity chromatography using Protein A. Single heavy and light chain bands on SDS-PAGE.
Biotinylation: Biotinylated and titrated for use in ELISA at 1/1000 dilution. Prepared in 1% BSA/50% glycerol/PBS, pH 7.4, 0.22µm filtered, preservative free.
Lot Number: xxxxx

Allergen Standard: Universal Allergen Standard
Composition: A formulation of eight purified natural allergens prepared in 1% BSA/50% glycerol/PBS, pH 7.4
Lot Number: xxxxx

Concentration:

Universal Allergen Standard	Protein Measurement	Concentration (ng/ml)
Der p 1	Amino-acid analysis	2500
Der f 1	Amino-acid analysis	2500
Der p 2	Amino-acid analysis	1000
Fel d 1	Amino-acid analysis	1000
Can f 1	Amino-acid analysis	2500
Rat n 1	Amino-acid analysis	1000
Mus m 1	Amino-acid analysis	250
Bla g 2	Amino-acid analysis	2500

ELISA Protocol for Mite Group 2 Allergens.

- Coat polystyrene microtiter plates (NUNC Maxisorp Cert. NUNC catalog # 439454) with 100µl mAb 1D8 at 10µl/10ml, i.e. 1/1000 dilution of stock, in 50mM carbonate-bicarbonate buffer, pH 9.6, incubate overnight at 4°C.
- Wash wells 3x with PBS-0.05% Tween 20, pH 7.4 (PBS-T). Incubate for 30 min. at room temperature with 100µl/well of 1% BSA, PBS-T. Wash 3x with PBS-T.
- Use doubling dilutions of the Universal Allergen Standard to make a control curve ranging from 100 - 0.2ng/ml Der p 2: Pipette 20µl Universal Allergen Standard into 180µl 1% BSA, PBS-T into wells A1 and B1 on the ELISA plate. Mix well and transfer 100µl across the plate into 100µl 1% BSA, PBS-T diluent to make 10 serial doubling dilutions. Wells A11, B11 and A12, B12 should contain only 1% BSA, PBS-T as blanks.
- Add 100µl of diluted allergen samples and incubate for 1 hour at room temperature. House dust extracts for Der p 2 analysis are routinely diluted two-fold from 1/10-1/80. Other sample types, like air filter extracts and allergen extracts, may require different dilutions.
- Wash wells 3x with PBS-T and add 100µl diluted biotinylated anti mAb 7A1. The antibody solution contains 50% glycerol and should be diluted 1/1000 in 1%BSA, PBS-T. Incubate for 1 hour at room temperature.
- Wash wells 3x and add 100µl diluted Streptavidin - Peroxidase (Sigma S5512, 0.25mg reconstituted in 1ml distilled water). The Streptavidin Peroxidase should be diluted 1/1000 in 1% BSA, PBS-T. Incubate for 30 minutes at room temperature.
- Wash wells 3x and develop the assays by adding 100µl 1mM ABTS in 70mM citrate phosphate buffer, pH 4.2 and 1/1000 dilution of H₂O₂. Read the plate when the absorbance at 405nm reaches 2.0-2.4.

Notes:

Universal Allergen Standard is recommended for immunoassay calibration purposes only. Not recommended for in-vitro antibody measurements, T cell studies, immunization purposes, or other uses.

Buffer recipes, storage conditions and a list of frequently asked questions can be found under "Protocols" on our web site: www.inbio.com.

For research and commercial use in vitro: not for human in vivo or therapeutic use.